## Vandprøvehenter

### Anvendelse

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## Water sampler

### Application

As the water sampler works according to the flow principle, the following must be observed during use:

1. Close both taps.
2. Let the water collector sink at its own weight vertically down to the desired sample depth. Never throw out the water collector as this may damage it.
3. To ensure there is only water from the sample depth in the water sampler, this is raised 3-5 times approx. 50 cm and quickly lowered again to the sample depth, after which the water sample collector is pulled up.
4. When draining, open the air screw - the round brass screw at the top of the cylinder - first.
5. If not already done: Fit the drain hose on the lower tap.
6. Hold the free end of the hose 1-3 mm above the bottom of the container to be used for analysis.
7. Open the drain tap (at the bottom of the cylinder).
8. For oxygen analysis, let the water sample fill the sample container from below and let 1-2 times the volume of the sample container float over. This ensures that the part of the water sample that was mixed with air during emptying through the hose is driven out of the sample holder.

NOTE: A sample container of 250 - 300 ml is best suited for the water sampler.

**Storage**

The water sampler should be stored horizontally or turned upside down. This allows it to dry completely, and the gaskets are not loaded. Avoid leaving the water sample collector lying in direct sunlight for extended periods of time. It can get very hot inside the container, and there is a risk of damaging the black rubber gaskets.

**Cleaning**

The water sample collector is cleaned by turning the bottom upside down and rinsing it with water.

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*© Frederiksen Scientific A/S*